


How to adjust the Face Shield

The visor is made in one piece and is put on easily with the neck part on the back of the head. Turn the round bubbles towards the head, but do not drag them to low on the forehead. The shield is made in one size, but the material is flexible and easy to adjust. To adjust the size of the shield, press the neckband on the sides or in the middle at the back of the neckband

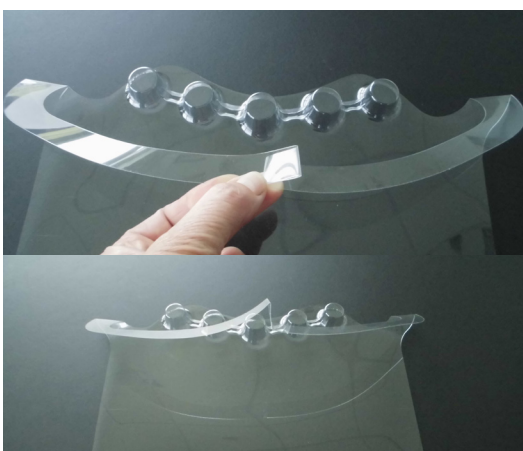

a) When bending the neck part towards you, in both sides before use, you create a self-adjustment spring-effect – leads to a one-size-fit function.



b) For XS/XXS head sizes you are able to bend one or two folds in center of the neck band, in order to adjust the neck band size.